

PureCube Ni-NTA Agarose

Cat. No.: 31103, 31105, 31110, 31112

Description	PureCube Ni-NTA Agarose is a resin for affinity purification of proteins fused to a polyhistidine tag (His-tag). A high density of nickel ions (Ni) loaded onto the ligand nitrilotriacetic acid (NTA) ensures a high binding capacity.
Support	7.5% agarose, highly cross-linked
Form	50% suspension in 20 mM sodium acetate with 20% ethanol, pH 6.6 (e.g. 10 ml will be 5 ml settled resin + 5 ml storage buffer)
Bead size	40 µm, spherical
Size exclusion limit	1 x 10 ⁶
Binding capacity	80 mg protein/ml resin. Binding capacity was determined with a 27 kDa 6xHis-tagged GFP protein. Please note that the binding capacity is protein dependent.
Maximum flow rate	≥1000 cm/h (at 15 cm bed height)
Maximum pressure	≥300 kPa (at 15 cm bed height)
Metal ion capacity	>15 µeqv Ni ²⁺ /ml
Compatible reagents	1 mM DTT, 1 mM EDTA, 100% methanol, 100% ethanol, 8 M urea, 6 M guanidinium hydrochloride, 30% (v/v) acetonitrile
Application	The suspension can be used for packing of gravity flow columns since the resin is optimized for column affinity chromatography, rather than for batch purification. To pack a column with 1 ml resin bed volume, 2 ml 50% suspension are necessary.
Stability	4 years after shipping
Storage	2-8 °C. Short-term: in neutral buffer (e.g. 50 mM phosphate, pH 7.0); Long-term: in neutral buffer with 20% ethanol
Shipping	Room temperature
Hazards	Product is not classified as hazardous according to (EC) No 1272/2008 [CLP]. A Material Safety Data Sheet is provided.
Manufacturer	Cube Biotech

For research use only

Trademark information

The owners of trademarks marked by "®" or "TM" are identified at <http://www.cube-biotech.com/information/patents>. Registered names, trademarks, etc. used in this document, even when not specifically marked as such, are not to be considered unprotected by law.

Important licensing information

This product is covered by intellectual property (IP) rights and on completion of the sale Cube Biotech grants respective Limited Use Label Licenses to purchaser. IP rights and Limited Use Label Licenses for said technology are further described and identified at <http://www.cube-biotech.com/information/patents> or upon inquiry at info@cube-biotech.com or at Cube Biotech GmbH, Creative-Campus-Allee 12, D-40789 Monheim, Germany. By use of this product the purchaser accepts the terms and conditions of all applicable Limited Use Label Licenses.